

### Quadriceps tendon tensile test

1. used wet cotton gauze to wrap the ends of sample and glued to both sides of clamps.
2. 1000 precondition cycles (except for SR protocol, image acquisition was every 10s, 0.1s for SR.)
3. 4.5 % max strain
4. 10%/s strain rate

oks00TR2-QAT-CXXX-09-01 lengths (mm): 12.4225 (for preconditioning), 12.6805 (for stress relaxation)

oks00TR2-QAT-CXXX-09-02 lengths (mm): 12.72 (for preconditioning), 12.859 (for stress relaxation)

oks00TR2-QAT-CXXX-09-03 lengths (mm): 12.853 (for preconditioning), 12.9655 (for stress relaxation)

